Pavella Creative **Design**

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BIOHUNGARICUM
From Genuine Source

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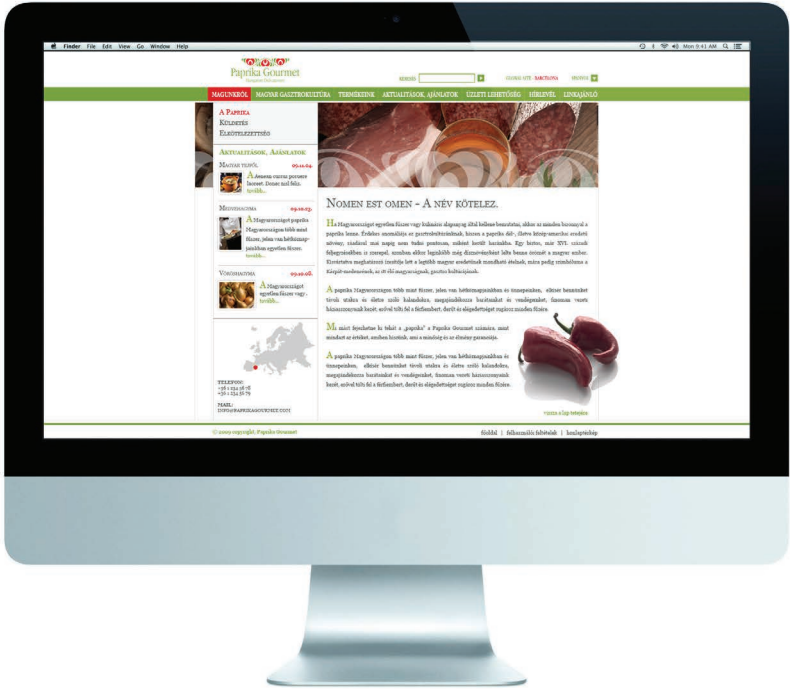
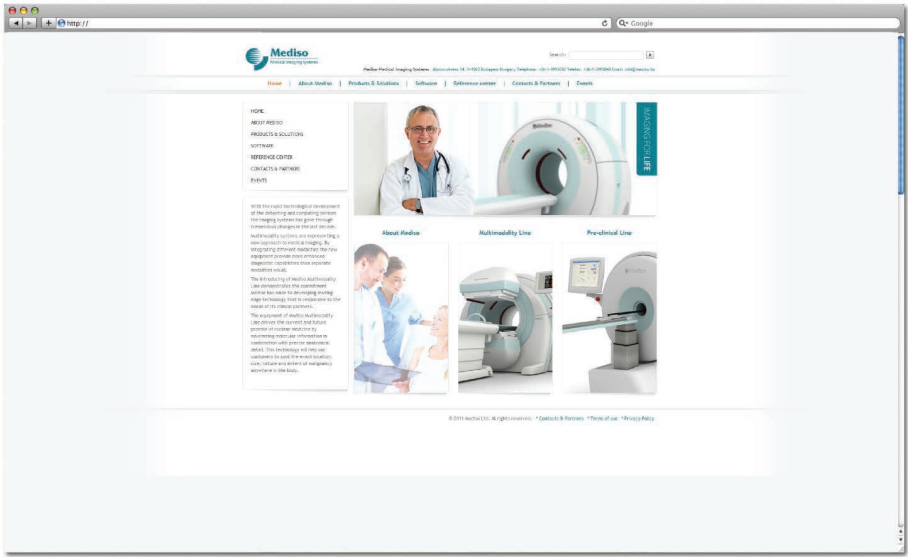
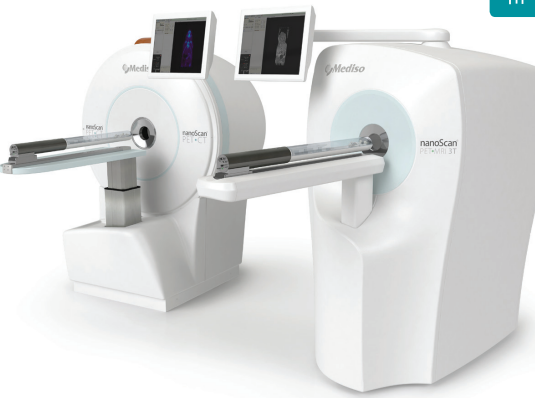
nanoScan® PET
Preclinical Imaging Systems

nanoScan® PET/MRI
nanoScan® PET/CT

IMAGING FOR SCIENCE

nanoScan® PC PET/CT
in vivo molecular and preclinical imager

The first sub-half mm³ PET volumetric resolution PET/CT with exceptional image quantification - accuracy over 97 %

Preclinical Line

IMAGING FOR SCIENCE

MEDISO GROUP OF COMPANIES

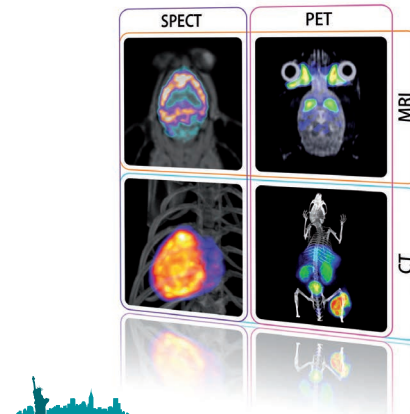
The most flexible clinical and preclinical imaging portfolio

IMAGING FOR LIFE

nanoScan® Family
the next generation four modality preclinical imaging platform
SPECT/MRI, PET/MRI, SPECT/CT, PET/CT, SPECT/CT/PET

nanoScan® - leading performance in all modalities

- breakthrough 13,000 cps/MBq SPECT sensitivity
- 300 µm SPECT and 700 µm PET resolution
- quantification accuracy over 97% for SPECT & PET
- Fully integrated SPECT/MRI and PET/MRI systems

IMAGING FOR LIFE

LEUCO-SCINT kit

Separation and labelling protocol (Aseptic conditions should be kept throughout the separation and labelling process)

1. Draw 2 ml of ACD solution and 3 ml of Fluorescein into each of the four 20 ml sterile plastic syringes.
2. Withdraw 15 ml patient's blood into each syringe and mix gently by inversion. Note in case of infants the labelling can be managed with 2-3 ml.
3. Leave to stand the syringes for 30-40 mins to allow red blood cells to sediment at room temperature.
4. When the red cells are sedimented to about half of the original volume of the blood, carefully draw up the leucocytes platelet rich plasma (LPP) into sterile tubes and centrifuge 10 mins at 150 g.
5. Remove and retain all the supernatant (platelet rich plasma, PRP). Pool the pellets of mixed leucocytes in one sterile tube. Centrifuge again for 10 mins at 150 g. Remove the remaining supernatant from the pellet of mixed leucocytes, leaving the pellet almost to dry. Note 1: During centrifugation phase start labelling a Leuco-Scint kit vial with 10-90s. Note 2: only 15 ml PRP is needed to procedure described in step 7.
6. Reconstitute one vial of LEUCO-SCINT with 1.5 ml eluate of ^{99m}Tc -generator containing 100-750 MBq ^{99m}Tc -pertechnetate. Shake for 10 mins to dissolve the HM-PAO and add exactly 1 ml immediately ^{99m}Tc -HM-PAO to a tube of mixed leucocytes (the activity of ^{99m}Tc -HM-PAO is approx. 400-500 MBq). Mix gently and incubate the cells for 10 mins at room temperature.
7. While incubating the leucocytes, centrifuge the PRP (step 5) for 5 mins at 2000 g to produce cell-free plasma (CFP).
8. Add 3.5 ml CFP (obtained in step 7) to the labelled cell suspension and gently mix. Centrifuge at 150g for 10 mins.
9. Transfer all supernatant to a tube.
10. Add 3.5 ml of CFP to the pellet of leucocytes, gently swirl to mix.
11. Measure the radioactivity and calculate the cell labelling efficiency (this activity of labelled leucocytes activity of labelled leucocytes ÷ activity of supernatant × 100).
12. The labelled leucocytes should be re-injected without delay. Imaging at 30 mins, 2 hrs and 24 hrs after administration.

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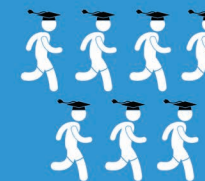
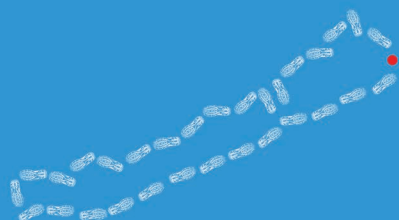
REFRESHMENTS ARE STILL
PRIMAVERA AND POWERBAR

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